




# ASEPTIQUIK® G SERIES DIMENSIONS

- A** = Height/Diameter
- B** = Total Length
- C** = Hose Barb Length

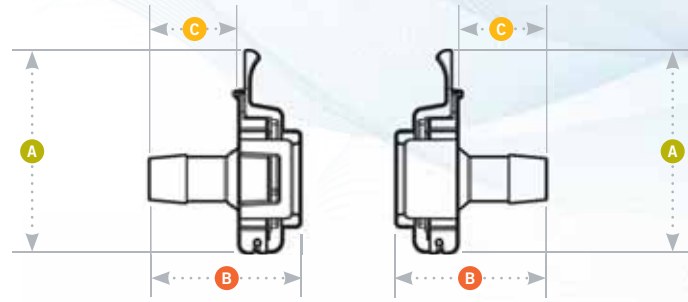

## Coupling Bodies

### • POLYCARBONATE with Blue Pull Tabs



#### TERMINATION

1/4" HOSE BARB  
3/8" HOSE BARB  
1/2" HOSE BARB  
3/4" HOSE BARB  
3/4" SANITARY

#### PART NO.

AQG17004  
AQG17006  
AQG17008  
AQG17012  
AQG33012

**A**

2.62  
2.62  
2.62  
2.62  
2.62

**B**

1.62  
1.62  
1.95  
2.36  
1.66

**C**

0.82  
0.82  
1.15  
1.56  
0.86

### • POLYCARBONATE HT with White Pull Tabs



#### TERMINATION

1/4" HOSE BARB  
3/8" HOSE BARB  
1/2" HOSE BARB  
3/4" HOSE BARB  
3/4" SANITARY

#### PART NO.

AQG17004HT  
AQG17006HT  
AQG17008HT  
AQG17012HT  
AQG33012HT

**A**

2.62  
2.62  
2.62  
2.62  
2.62

**B**

1.62  
1.62  
1.95  
2.36  
1.66

**C**


0.82  
0.82  
1.15  
1.56  
0.86



4040 Norex Drive, Chaska MN 55318  
phone: 952.448.4412 | fax: 952.448.6180  
sales@smithengineeringinc.com  
www.smithengineeringinc.com


Validation reports  
are available at  
[cpcworldwide.com/bio](http://cpcworldwide.com/bio).

## FLIP-CLICK-PULL Assembly Procedure

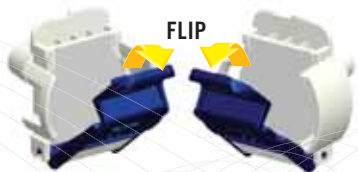


**1**

Unsnap and flip down the protective pull tab covers on each connector half.

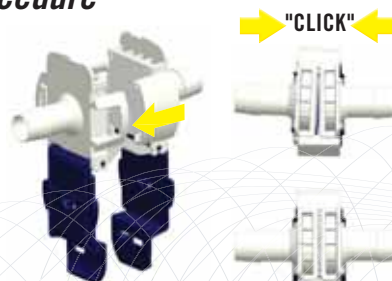

**2**

Align the connector halves with the pull tabs hanging down. Slide the two halves together. Independently squeeze each side of the connector until you hear an audible "CPC Click."

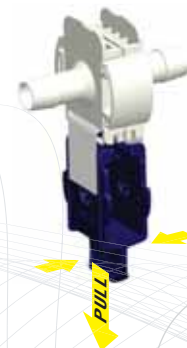

**3**

To complete the connection, simply snap the pull tabs together by pushing on the CPC logos and pull the membranes from the connector.